

## Deodorant

(Powder)

Sample Category: Miscellaneous

### Sample Requirements

Sample Weight: 0.2g

Reagents: 2mL Conc. HNO<sub>3</sub>, 6mL HCl

Vessels: 35mL glass, quartz, and Teflon™ lined

### Microwave Heating Program

Control Style: Ramp To Temperature




Pressure Mode: Inorganic

### Recommended Equipment



Discover SP-D

Stage	Temp (°C)	Ramp Time	Hold Time	Pressure (psi)	Power (W)	Stirring
1	210	5:00	5:00	400	300	Med

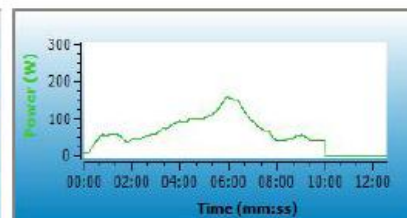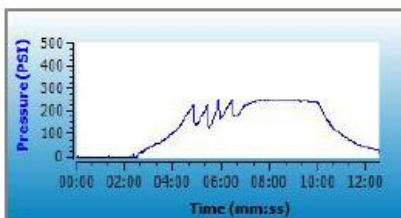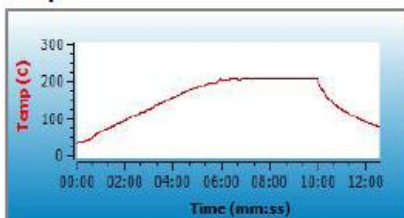

#### Notes:

1. If using a vial other than the recommended choice, adjust sample size and pressure limit to values appropriate for the vessel chosen.

#### General Precautions:

This procedure is a reference for sample digestion using a CEM System and may need to be modified or changed to obtain the required results on your sample