

**PVC**  
(film)

Sample Category: Oils and Plastics

**Sample Requirements**

Sample Weight: 0.2g  
 Reagents: 10mL Conc. HNO<sub>3</sub>  
 Vessels: 35mL glass, quartz, and Teflon™ lined

**Microwave Heating Program**



Control Style: Ramp To Temperature  
 Pressure Mode: Organic

**Recommended Equipment**



Discover SP-D

Stage	Temp (°C)	Ramp Time	Hold Time	Pressure (psi)	Power (W)	Stirring
1	210	5:00	3:00	400	300	Med

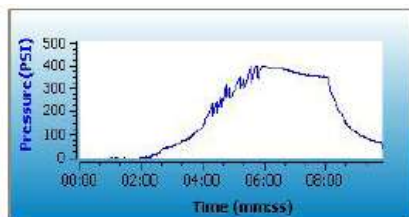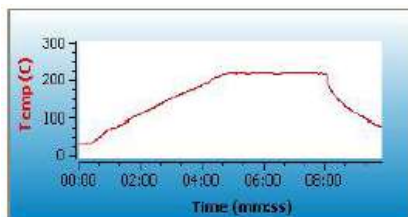

**Notes:**

1. If using a vial other than the recommended choice, adjust sample size and pressure limit to values appropriate for the vessel chosen.

**General Precautions:**

This procedure is a reference for sample digestion using a CEM System and may need to be modified or changed to obtain the required results on your sample