

## Sulfuric Char

Sample Category: Oils and Plastics

### Sample Requirements

Sample Weight: 0.2g  
 Reagents: 6mL Conc. H<sub>2</sub>SO<sub>4</sub>  
 Vessels: 35mL glass, quartz, and Teflon™ lined

### Recommended Equipment






Discover SP-D

### Microwave Heating Program

Control Style: Ramp To Temperature  
 Pressure Mode: inorganic

Stage	Temp (°C)	Ramp Time	Hold Time	Pressure (psi)	Power (W)	Stirring
1	240	6:00	3:00	400	300	Med

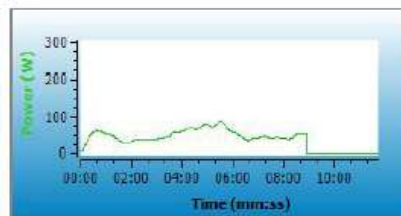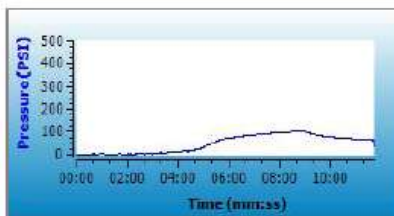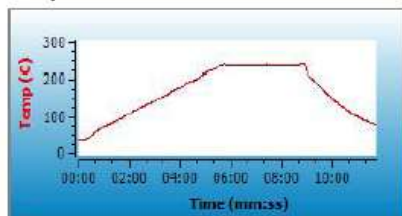

### Notes:

1. If using a vial other than the recommended choice, adjust sample size and pressure limit to values appropriate for the vessel chosen.

### General Precautions:

This procedure is a reference for sample digestion using a CEM System and may need to be modified or changed to obtain the required results on your sample