

## Tomato Leaves

Sample Category: Food / Plant / Animal Tissue

### Sample Requirements

Sample Weight: 0.3g

Reagents: 5mL Conc. HNO<sub>3</sub>, 2mL Deionized Water, 1mL Conc. HF

Vessels: 35mL glass, quartz, and Teflon™ lined

### Recommended Equipment





Discover SP-D

### Microwave Heating Program

Control Style: Ramp To Temperature

Pressure Mode: Organic

Stage	Temp (°C)	Ramp Time	Hold Time	Pressure (psi)	Power (W)	Stirring
1	200	4:00	2:00	400	300	Med

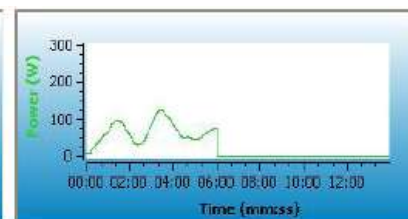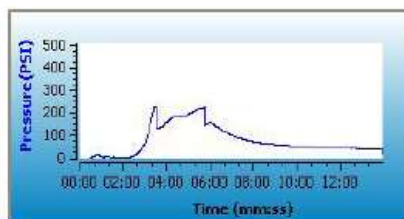

#### Notes:

1. If using a vial other than the recommended choice, adjust sample size and pressure limit to values appropriate for the vessel chosen.
2. The addition of Deionized water may improve solubility of minerals and prevent temperature spikes due to exothermic reactions.

#### Safety Precautions for Working with Samples Requiring HF

1. The instrument must be placed in a hood and the SP-D HF Scrubber must be installed.
2. Sample mass must be  $\leq 0.3$  grams for organic samples and  $\leq 0.5$  grams for inorganic samples.
3. The volume of HF required for total digestion may vary depending upon the sample matrix. Do not exceed 3 mL of total HF volume in the 35 mL Teflon lined vials.
4. It is recommended that Personal Protective Equipment (PPE) for handling HF include chemical splash goggles with a safety shield, lab coat with chemical splash apron, and heavy weight nitrile or rubber gloves.

#### General Precautions:

This procedure is a reference for sample digestion using a CEM System and may need to be modified or changed to obtain the required results on your sample

