



Procedure

Weigh 0.1 g of sample into the digestion vessel. Add 3 mL of H₃PO₄, 2 mL of HNO₃, and 3 mL of HBF₄. HBF₄ should be added slowly and carefully to the sample. Allow any initial reaction to subside before sealing the vessel.

Notes

This application can only be run in the iPrep vessel.

Recommended Equipment

MARS 6 iWave

Recommended Vessels

iPrep

Reagents

H₃PO₄
HNO₃
HBF₄

Max Sample Weight

0.1 g

Sample Type

Organic

Control Type

Ramp to Temperature

Method Type

One Touch

Heating Program

Stage	Temp (°C)	*Ramp (mm:ss)	Hold (mm:ss)	Pressure (psi)	* Power (W)	Stirring
1	230	30:00	30:00	N/A	700-1800	Off

* Ramp times and power may vary depending on the type and number of vessels.

Results

Sample was clear, colorless, and particle free upon dilution to 50 mL.

General Precaution

- This procedure is a reference point for sample digestion using a CEM system and may need to be modified or changed to obtain the required results on your sample.
- If using a vessel other than the recommended choice, adjust sample size and pressure limit to values appropriate for the vessel chosen.
- Manual venting of CEM vessels should be performed when wearing hand/eye/body protection and when the vessel contents are at or below room temperature to avoid the potential for chemical burns. Always point the vent hole away from the operator.
- If programming as One Touch, the ramp time and power will be automatically determined based on the number and type of vessels detected.