



Systems

SMART 6

Equipment

CEM Square Glass Fiber Pads
Disposable Pipette
Sample Press

SMART 6 Settings

| Calculation Mode | Option | Display Mode |
|------------------|----------|--------------|
| Moisture/Solids | Standard | Solids |

SMART 6 Categories

| Ramp Time (sec) | Delta Weight (mg) | Delta Time (sec) |
|-----------------|-------------------|------------------|
| 30 | 5 | 20 |

| Am. Fan Speed (%) | Run Fan Speed (%) | Run Fan Off (sec) | Weight Comp. (On /Off) |
|-------------------|-------------------|-------------------|------------------------|
| 0 | 50 | 60 | On |

Parameters

| Stage | Mag (%) | IR (%) | Time (mm:ss) | Max Temp (°C) | Mode |
|-------|---------|--------|--------------|---------------|--------------|
| 1 | 100 | 50 | 10:00 | 110 | Constant Wt. |

Sample Preparation

Procedure

Spread Technique