



#### Systems

SMART Q

#### Equipment

CEM Square Glass Fiber Pads  
Disposable Pipette

#### SMART 6 Settings

Calculation Mode	Option	Display Mode
Moisture/Solids	Standard	Moisture

#### SMART 6 Categories

Ramp Time (sec)	Delta Weight (mg)	Delta Time (sec)
0	5	20

Am. Fan Speed (%)	Run Fan Speed (%)	Run Fan Off (sec)	Weight Comp. (On /Off)
0	50	30	On

#### Parameters

Stage	IR (%)	Time (mm:ss)	Max Temp (°C)	Mode
1	100	10:00	100	Constant Wt.

#### Sample Preparation

Mix sample thoroughly.

#### Procedure

Tare two glass fiber sample pads.  
Pipette 1.9-2.1 grams of sample on one pad and cover with the second pad.  
Place the pads with sample centered on balance pan. Close cover, and press "START".

#### Spread Technique