



### Systems

SMART 6

### Equipment

CEM Square Glass Fiber Pads  
Disposable Pipette  
Sample Press

### SMART 6 Settings

Calculation Mode	Option	Display Mode
Moisture/Solids	Standard	Solids

### SMART 6 Categories

Ramp Time (sec)	Delta Weight (mg)	Delta Time (sec)
60	2	20

Am. Fan Speed (%)	Run Fan Speed (%)	Run Fan Off (sec)	Weight Comp. (On /Off)
0	100	30	On

### Parameters

Stage	Mag (%)	IR (%)	Time (mm:ss)	Max Temp (°C)	Mode
1	100	100	10:00	150	Constant Wt.

### Sample Preparation

Mix sample thoroughly.

### Procedure

Tare two glass fiber sample pads. Apply 3.9-4.1 grams of sample on one pad and cover with the second pad. Press pads containing sample with sample press. Place the pads with sample centered on balance pan. Close cover, and press "START".

### Spread Technique