

# Routine Maintenance Log

## CEM SMART Turbo & SMART Trac II



		Daily	Weekly		Monthly		PM
		SMART Trac II	SMART Turbo	SMART Trac II	SMART Turbo	SMART Trac II	
		Perform a Frequency Optimization	Clean the SMART Turbo interior Inspect the SMART Turbo Air Shield Verify Operation of Energy Distributor Blade	Perform a Pulse Width Optimization	Perform Balance Calibration Perform MONITOR Sample Test Clean the Air Intake Vents	Perform a Check Standard Inspect the Air filters	Performed by CEM Service
J F M A M J J A S O N D	1	<input type="checkbox"/>					
	2	<input type="checkbox"/>					
	3	<input type="checkbox"/>					
	4	<input type="checkbox"/>					
	5	<input type="checkbox"/>					
	6	<input type="checkbox"/>					
	7	<input type="checkbox"/>	<input type="checkbox"/>	<input type="checkbox"/>	<input type="checkbox"/>		
	8	<input type="checkbox"/>					
	9	<input type="checkbox"/>					
	10	<input type="checkbox"/>					
	11	<input type="checkbox"/>					Date
	12	<input type="checkbox"/>					_____
	13	<input type="checkbox"/>					
	14	<input type="checkbox"/>	<input type="checkbox"/>	<input type="checkbox"/>	<input type="checkbox"/>		
	15	<input type="checkbox"/>					
	16	<input type="checkbox"/>					CEM Technician
	17	<input type="checkbox"/>					_____
	18	<input type="checkbox"/>					
	19	<input type="checkbox"/>					
	20	<input type="checkbox"/>					
	21	<input type="checkbox"/>	<input type="checkbox"/>	<input type="checkbox"/>	<input type="checkbox"/>		
	22	<input type="checkbox"/>					
	23	<input type="checkbox"/>					
	24	<input type="checkbox"/>					
	25	<input type="checkbox"/>					
	26	<input type="checkbox"/>					
	27	<input type="checkbox"/>					
	28	<input type="checkbox"/>	<input type="checkbox"/>	<input type="checkbox"/>	<input type="checkbox"/>	<input type="checkbox"/>	<input type="checkbox"/>
	29	<input type="checkbox"/>					
	30	<input type="checkbox"/>					
	31	<input type="checkbox"/>					

**Daily**

- Perform a frequency optimization

**Weekly**

- Clean the SMART Turbo interior
- Inspect the air shield and replace if necessary
- Verify proper operation of the energy distributor blade
- Perform a pulse width optimization

**Monthly**

- Perform a balance calibration
- MONITOR sample test
- SMART Turbo air intake vents
- Perform a check standard
- Inspect the air filters and clean if necessary

**Tool Requirements**

- 10 gram weight
- Glass cleaner
- Lint-free towels
- Small soft bristle brush

**CEM Contacts**

**Service**  
Email: [service@cemhelp.com](mailto:service@cemhelp.com)  
Phone: 800-726-5551

**Sales**  
Email: [sales@cem.com](mailto:sales@cem.com)  
Phone: 800-726-3331

**Laboratories**  
Email: [labs@cem.com](mailto:labs@cem.com)  
Phone: 800-726-3331

CEM Corporation  
3100 Smith Farm Road  
Matthews, NC 28104

Refer to the routine maintenance instructions outlined in the SMART Turbo and SMART Trac II Operation Manuals for a more comprehensive explanation of the requirements. This document is intended as a daily reference and record of required routine maintenance activity.

### Preventive Maintenance (PM)

Preventive maintenance is performed by a CEM Certified Technician, typically once per year. However, should individual requirements vary, a CEM representative is available to assist in creating a schedule to meet your maintenance needs. Contact the CEM Contract Administrator at (800) 726-3331 for assistance.

