

Moisture Analysis of Milk

M601.0

Parameters

Category: Dairy

Stage 1: Constant Weight

iPower Mag: 100%

iPower IR: 000%

Temperature: 100°C

Weight Compensation: ON

Recommended Systems

SMART 6

Equipment

CEM Square Glass Fiber Sample Pads

Pipette

Sample Preparation

Stir well before analysis.

Procedure

Tare 2 glass fiber sample pads. With the pads still on the balance pan, puddle 3 grams of sample in the middle of the top pad, do not cover. Place the pads containing the sample on the balance pan, close cover and press START.

Precision Range:

Moisture $\pm 0.05\%$

*The above parameters are suggested starting points and may need to be modified to meet certain product specifications