

Digestion of Yellow Pill (whole)

Sample Category: Pharmaceuticals

Recommended Equipment

Sample Requirements

Sample Weight: 1 g

Reagents: 10 mL Conc. HNO₃

2 mL H₂O₂ (30%)

Vessels: 80 mL quartz.

Microwave Heating Program

Note: Add acid slowly and allow sample to pre-digest at Least 10 minutes before running




Control Style: Ramp To Temperature

Pressure Mode: Organic



Discover SP-D 80 ml

Stage	Temp (°C)	Ramp Time	Hold Time	Pressure (psi)	Power (W)	Stirring
1	200	05:00	03:00	400	300	Med

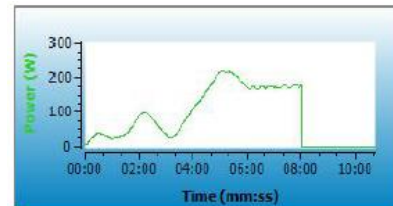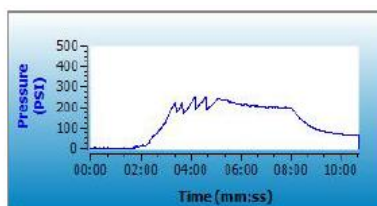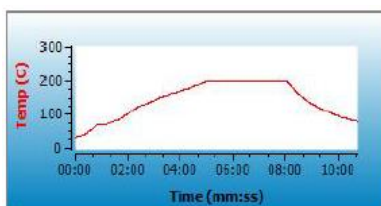

Notes:

1. If using a vial other than the recommended choice, adjust sample size and pressure limit to values appropriate for the vessel chosen.
2. This method may not provide a total digest of all sediment samples.

General Precautions:

This procedure is a reference for sample digestion using a CEM System and may need to be modified or changed to obtain the required results on your sample