

EDGE Release Notes


Software v1.23

In order to update to v1.23, v1.22 must first be installed.

Software Features Added

- Volume accuracy for low volatile solvents improved
- Power Limiting Temperature

This feature can be found by selecting System Menu -> Settings -> Run and the recommended value is 71°C. This feature provides better heater control for temperatures below the entered value and reduces temperature overshoot errors. This feature is only for a method temperature of 100°C or below.

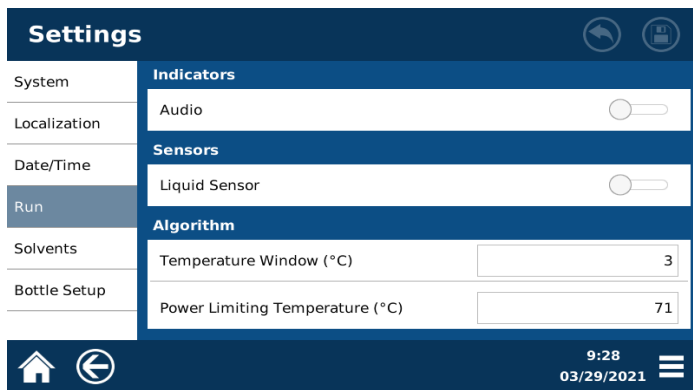

Software v1.22

Q-Disc Updates

S1 (Part# 153530): New packaging with preassembled G1 Q-Disc sandwiched between two C9 Q-Discs

If currently using G1, order these when you run out of G1 and be aware that they contain 125 Q-Discs in a box instead of 250. You will have to change the Q-Disc in your method to S1 which will result in losing run data. If this is a concern, copy this method and select S1 as the Q-Disc. As long as you do not edit any of the parameters from the original method, your run data will be saved.

