



Procedure

Weigh 0.1 g of the sample into the digestion vessel. Add 8 mL H₃PO₄. Gently swirl the mixture and wait approximately 15 minutes to predigest before closing the vessel.

Notes

This application can only be run in the iPrep vessel.

This method may not provide a total digest of all samples. Hydrofluoric acid will be required to provide complete digestion of some sample matrixes.

Recommended Equipment

MARS 6 iWave

Recommended Vessels

iPrep

Reagents

H₃PO₄

Max Sample Weight

0.5 g

Sample Type

Organic

Control Type

Ramp to Temperature

Method Type

One Touch

Heating Program

Stage	Temp (°C)	*Ramp (mm:ss)	Hold (mm:ss)	Pressure (psi)	* Power (W)	Stirring
1	260	25:00	45:00	N/A	1800	Off

* Ramp times and power may vary depending on the type and number of vessels.

Results

Sample was clear, colorless with some white particles remaining upon dilution to 50 mL.

General Precaution

- This procedure is a reference point for sample digestion using a CEM system and may need to be modified or changed to obtain the required results on your sample.
- If using a vessel other than the recommended choice, adjust sample size and pressure limit to values appropriate for the vessel chosen.
- Manual venting of CEM vessels should be performed when wearing hand/eye/body protection and when the vessel contents are at or below room temperature to avoid the potential for chemical burns. Always point the vent hole away from the operator.
- If programming as One Touch, the ramp time and power will be automatically determined based on the number and type of vessels detected.