



Systems

ORACLE
SMART 6

Equipment

CEM Square Glass Fiber Pads
CEM Trac Film
ORACLE Station
Sample Press
Spatula

SMART 6 Settings

Calculation Mode	Option	Display Mode
Moisture/Fat	Standard	Moisture

SMART 6 Categories

Ramp Time (sec)	Delta Weight (mg)	Delta Time (sec)
60	2	20

Am. Fan Speed (%)	Run Fan Speed (%)	Run Fan Off (sec)	Weight Comp. (On /Off)
0	50	30	Off

ORACLE Parameters

Cooling Enabled	Temperature (°C)	QuickPrep™ Temp (°C)	QuickPrep™ Time (sec)	Warm Delay (sec)	Max Cool Time
Yes	45	70	15	5	3:30

Parameters

Stage	Mag (%)	IR (%)	Time (mm:ss)	Max Temp (°C)	Mode
1	100	0	3:30	140	Set Time

Sample Preparation

Blend sample using an industrial or consumer grade food processor until homogeneous.

Procedure

Spread Technique

