



**Systems**

SMART Q

**Equipment**

CEM Pad Press  
CEM Square Glass Fiber Pads  
Spatula

**SMART 6 Settings**

Calculation Mode	Option	Display Mode
Moisture/Solids	Standard	Moisture

**SMART 6 Categories**

Ramp Time (sec)	Delta Weight (mg)	Delta Time (sec)
0	2	20

Am. Fan Speed (%)	Run Fan Speed (%)	Run Fan Off (sec)	Weight Comp. (On /Off)
0	50	30	On

**Parameters**

Stage	IR (%)	Time (mm:ss)	Max Temp (°C)	Mode
1	100	8:30	150	Set Time

**Sample Preparation**

Mix sample thoroughly.

**Procedure**

Prepare two glass fiber sample pads with the CEM Pad Press and tare. Spread 10-20 grams of sample across the pad indentation with a spatula. Place the pad with sample centered on balance pan. Close cover, and press "START".

**Spread Technique**