



Systems

SMART Q

Equipment

CEM Basket
CEM Square Glass Fiber Pads
Spatula

SMART 6 Settings

Calculation Mode	Option	Display Mode
Moisture/Solids	Standard	Moisture

SMART 6 Categories

Ramp Time (sec)	Delta Weight (mg)	Delta Time (sec)
0	2	20

Am. Fan Speed (%)	Run Fan Speed (%)	Run Fan Off (sec)	Weight Comp. (On /Off)
0	50	30	On

Parameters

Stage	IR (%)	Time (mm:ss)	Max Temp (°C)	Mode
1	100	10:00	180	Set Time

Sample Preparation

None.

Procedure

Use CEM Basket lined with one glass fiber pad. Tare lined basket, then spread 14.9-15.1 grams of sample into basket, place on the balance pan. Close cover, and press "START

Spread Technique